



Emergency Eyewash/Drench Hose Inspection Checklist

Building: _____ Room #: _____

How to Test:

- Visually inspect the area around the eyewash/drench hose to ensure the area is clear of obstructions and all protective covers are in good condition;
- Activate and flush eyewash/drench hose for 3 minutes or until water is clear ;
- Ensure all jets are working properly with good flow;
- Ensure the eyewash/drench hose water is clear before turning off;
- If there is an issue with your eyewash you can contact facilities or EH&S for assistance; and
- Refer to the [Eyewash & Safety Shower Safety Instruction](http://oregonstate.edu/ehs/sites/default/files/pdf/si/eyewash_and_safety_shower_si.pdf) for more information:
http://oregonstate.edu/ehs/sites/default/files/pdf/si/eyewash_and_safety_shower_si.pdf

Date	Initials

Date	Initials

Date	Initials

Questions/Concerns: Contact EH&S
Office Phone: 541-737-2273
Website: ehs.oregonstate.edu